



cGK Positive Control (catalytic domain)

Product Data Sheet

For Research Use Only, Not for use in diagnostic procedures

## cGMP Dependent Protein Kinase Positive Control (catalytic domain)

(Human, a.a. 324-671, recombinant protein expressed in Sf9)

Cat# CY-E1161-1

Lot No.  
For 200 Assays  
(40 units /  $\mu\text{L}$  x 100 $\mu\text{L}$ )

**Product Description:** Human catalytic domain of cGMP-Dependent Protein Kinase containing a N-terminal GST tag and a C-terminal His tag, expressed in recombinant baculovirus infected sf9 cells. Purified by sequentially using GSH agarose and Ni-NTA agarose chromatography. The GMP Dependent Protein Kinase Positive Control (catalytic domain) is designed to use for CycLex cGK Assay Kit [Cat# CY-1161]. The GMP Dependent Protein Kinase Positive Control (catalytic domain) should be added to the well at 2 units/well. For instance, diluted positive control 1:200, use 10  $\mu\text{L}$  for 1 assay. Unused GMP Dependent Protein Kinase Positive should be stored at  $-70^{\circ}\text{C}$ .

The GMP Dependent Protein Kinase Positive Control (full length) [Cat# CY-E1161-2] is also available from CycLex.

**Product Size:** Catalytic domain of cGMP-Dependent Protein Kinase (I alpha-Isozyme) 4000 units/100  $\mu\text{L}$

**Formulation:** The GMP Dependent Protein Kinase Positive Control is supplied frozen in a buffer containing 20mM Hepes-KOH (pH 7.5), 1 % BSA, 1mM EDTA, 1 mM DTT, 50mM NaCl, 0.03 % Brij35 and 50% glycerol.

**Source:** Recombinant cGMP-Dependent Protein Kinase catalytic domain is purified from recombinant baculovirus infected Sf9 cells. This enzyme shows protein kinase activity in the absence of cGMP under the protocol described below.

**Molecular Weight:** cGMP-Dependent Protein Kinase demonstrates a single 64 kDa band by SDS-PAGE analysis.

**Purity:** cGMP-Dependent Protein Kinase is greater than 90% pure as determined by SDS-PAGE analysis.

**Substrates:** cGMP-Dependent Protein Kinase phosphorylates a number of substrates, including histone proteins H2b and H4, brain G protein and high mobility group 14 protein.

**Inhibitors:** Polycations, such as poly-L-arginine, inhibit cGMP-Dependent Protein Kinase.

**Unit Definition:** One unit is defined as the amount of kinase required to incorporate 1pmol of phosphate into the GST-G substrate fusion protein, per minute at  $30^{\circ}\text{C}$ .

**Assay Conditions:** Assay activity of cGMP-Dependent Protein Kinase (catalytic domain) in a 50  $\mu\text{L}$  reaction containing 20 mM Hepes KOH (pH 7.5), 5 mM  $\text{MgCl}_2$ , 1 mM DTT, 100  $\mu\text{M}$  [ $\gamma$ - $^{32}\text{P}$ ] ATP (1  $\mu\text{Ci}$ ) and 4  $\mu\text{g}$  of GST-G substrate fusion protein in the absence of cGMP. Start the reaction by adding



cGK Positive Control (catalytic domain)

Product Data Sheet

**For Research Use Only, Not for use in diagnostic procedures**

10 $\mu$ L of the enzyme, diluted 50-fold in a buffer containing 20 mM Hepes KOH (pH 7.5), 1 mM DTT, 0.03 % Brij35. Incubate for 30 minutes at 30°C. Terminate the reaction by adding 600  $\mu$ L of cold 10 % TCA solution containing 0.2 % sodium pyrophosphate and stand on ice for 15 min. Filterate acid insoluble material through GFC filters (Whatman Inc.), wash 4 times with 1 % TCA and rinse filters with ethanol. Dry filters and count in a liquid scintillation counter.

**Storage Conditions:** Store cGMP-Dependent Protein Kinase frozen at  $-70^{\circ}\text{C}$ , where it is stable for at least 12 months from date of purchase. Avoid multiple freeze-thaw cycles. If thawed, store at  $4^{\circ}\text{C}$  if entire vial will be used within two weeks; or refreeze aliquots in liquid nitrogen and store at  $-70^{\circ}\text{C}$ .

**Related Products:**

\*Cyclic GMP dependent protein kinase (cGK Assay Kit): Cat# CY-1161

\*cGMP Dependent Protein Kinase Positive Control (full length): Cat# CY-E1161-2

**References**

1. Edelman, A.M., Blumenthal, D.K. and Krebs, E.G. Protein serine/threonine kinases. *Ann. Rev. Biochem.* **56**, 567, 1987
2. Beebe, S.J. and Corbin, J.D. In: *The Enzymes*, Vol. **17**, 3rd ed., Boyer, P.D. and Krebs, E.D., eds., 44, 1986
3. Corbin, J.D. and Doskeland, S.O. *J. Biol. Chem.* **258**, 11391-1983
4. Endo, S et al. *Proc. Natl. Acad. Sci. USA.* **96**, 2467-2472, 1999

**PRODUCED BY**

CycLex Co., Ltd.

1063-103 Terasawaoka

Ina, Nagano 396-0002

Japan

Fax: +81-265-76-7618

e-mail: [info@cyclex.co.jp](mailto:info@cyclex.co.jp)

URL: <http://www.cyclex.co.jp>

CycLex/CircuLex products are supplied for research use only. CycLex/CircuLex products and components thereof may not be resold, modified for resale, or used to manufacture commercial products without prior written approval from CycLex Co., Ltd. To inquire about licensing for such commercial use, please contact us via email.