



Plk1 Positive Control

Product Data Sheet

For Research Use Only, Not for use in diagnostic procedures

## Plk1 Positive Control

(Human, full length, recombinant enzyme expressed in *E. coli*.)

Cat# CY-E1163

Lot No.

For 200 Assays

(0.02 units /  $\mu\text{L}$  x 100  $\mu\text{L}$ )

**Product Description:** Constitutive active form of human full length Plk1, in which there is one mutation T210D in an active loop of kinase domain, containing a *N*-terminal GST tag, expressed in *E. coli*. Purified by GSH agarose chromatography. The Plk1 Positive control is designed to use for Plk1 Assay/Inhibitor Screening Kit [Cat# CY-1163]. The Plk1 Positive Control should be added to the well at 0.1 m units/well. For instance, diluted positive control 1:2000, use 10  $\mu\text{L}$  for 1 assay. Unused Plk1 Positive control should be stored at  $-70^{\circ}\text{C}$ .

**This product is produced under license: U.S. Pat. No. 6,180,380**

**Product Size:** Full length Plk1: 2 units/100  $\mu\text{L}$

**Formulation:** The Plk1 Positive Control is supplied frozen in a buffer containing 20mM Hepes-KOH (pH 7.5), 1 % BSA, 1mM EDTA, 1 mM DTT, 50mM NaCl, 0.03 % Brij35 and 50% glycerol.

**Source:** Human full length Plk1, containing *N*-terminal GST tag, expressed in *E. coli*.

**Molecular Weight:** Plk1 demonstrates a single 92 kDa band by SDS-PAGE analysis.

**Purity:** Plk1 is greater than 85 % pure as determined by SDS-PAGE analysis.

**Substrates:** Plk1 phosphorylates a number of substrates, including Cyclin B, MBP.

**Inhibitors:** Specific Plk1 inhibitor has not been discovered yet.

**Unit Definitions:** One unit is defined as the amount of kinase required to incorporate 1 nmol of phosphate into alpha-casein per minute at  $30^{\circ}\text{C}$ .

**Assay Conditions:** Assay activity of Plk1 in a 50  $\mu\text{L}$  reaction containing 20 mM Hepes KOH (pH 7.5), 5 mM  $\text{MgCl}_2$ , 1 mM DTT, 100  $\mu\text{M}$  [ $\gamma$ - $^{32}\text{P}$ ] ATP (1  $\mu\text{Ci}$ ), and 4  $\mu\text{g}$  of alpha-casein. Start the reaction by adding 10 $\mu\text{L}$  of the enzyme, diluted 50-fold in a buffer containing 20 mM Hepes KOH (pH 7.5), 1 mM DTT, 0.03 % Brij35. Incubate for 30 minutes at  $30^{\circ}\text{C}$ . Terminate the reaction by adding 600  $\mu\text{L}$  of cold 10 % TCA solution containing 0.2 % sodium pyrophosphate and stand on ice for 15 min. Filtrate acid insoluble material through GFC filters (Whatman Inc.), wash 4 times with 1 % TCA and rinse filters with ethanol. Dry filters and count in a liquid scintillation counter.

**Storage and Stability:** Stable for 12 months at  $-70^{\circ}\text{C}$  from date of shipment. For maximum recovery of product, centrifuge the original vial after thawing and prior to removing the cap. Aliquot enzyme to avoid repeated freezing and thawing.

### Related Products:

\*Plk1 Assay/Inhibitor Screening Kit: Cat# CY-1163



Plk1 Positive Control

Product Data Sheet

**For Research Use Only, Not for use in diagnostic procedures**

**General References:**

1. Lane, H. A. & Nigg, E. A. *Trends Cell Biol.* **7**, 63-68, 1997
2. Glover, D. M., Hagan, I. M. & Tavares, A. A. M. *Genes Dev.* **12**, 3777-3787, 1998
3. Sunkel, C. E. & Glover, D. M. *J. Cell Sci.* **89**, 25-38, 1988
4. Ohkura, H., Hagan, I. & Glover, D. M. *Genes Dev.* **9**, 1059-1073, 1995
5. Song, S. & Lee, K. S. *J. Cell Biol.* **152**, 451-469, 2001
6. Lane, H. A. & Nigg, E. A. *J. Cell Biol.* **135**, 1701-1713, 1996
7. Kotani, S., Tugendreich, S., Fujii, M., Jorgensen, P., Watanabe, N., Hoog, C., Hieter, P. & Todokoro, K. *Mol. Cell* **1**, 371-380, 1998
8. Alexandru, G., Uhlmann, F., Mechtler, K., Poupart, M. & Nasmyth, K. *Cell* **105**, 459-472, 2001
9. Liu, X.; Erikson, R. L: Polo-like kinase (Plk1) depletion induces apoptosis in cancer cells. *Proc. Nat. Acad. Sci.* **100**: 5789-5794, 2003.

**This product is produced under license: U.S. Pat. No. 6,180,380**

**PRODUCED BY**

CycLex Co., Ltd.  
1063-103 Terasawaoka  
Ina, Nagano 396-0002  
Japan  
Fax: +81-265-76-7618  
e-mail: [info@cyclex.co.jp](mailto:info@cyclex.co.jp)  
URL: <http://www.cyclex.co.jp>

CycLex/CircuLex products are supplied for research use only. CycLex/CircuLex products and components thereof may not be resold, modified for resale, or used to manufacture commercial products without prior written approval from CycLex Co., Ltd.. To inquire about licensing for such commercial use, please contact us via email.